

Technical Specifications

EIGER2 S 9M

Document Version v1.6.1

DECTRIS Ltd.

5405 Baden-Daettwil
Switzerland
www.dectris.com

CONTENT

CONTENT	i
DOCUMENT HISTORY	ii
Current Document	ii
Changes	ii
1 GENERAL INFORMATION	1
1.1 Contact and Support	1
1.2 Explanation of Symbols	1
1.3 Warranty Information	2
1.4 Disclaimer	2
2 USE OF THE EIGER2 S 9M	3
2.1 Product Return and Recycling	3
3 TECHNICAL SPECIFICATIONS	4
3.1 Specifications	4
3.1.1 Detector	4
3.2 Ratings	5
3.2.1 Detector	5
3.2.2 Detector Control Unit	6
3.2.3 Thermal Stabilization Unit	6
3.3 Ambient Conditions	7
4 DETECTOR DIMENSIONS AND CONNECTORS	8
4.1 EIGER2 S 9M Detector	8
4.1.1 Technical Drawing	8
4.1.2 Front Side of the Detector	9
4.1.3 Back Side of the Detector	10
4.1.4 Status LEDs	11
4.1.5 Connectors and Connecting Cables/Pipes	11
4.2 Detector Control Unit	12
4.2.1 Configuration	12
4.2.2 Connectors	14
4.3 Thermal Stabilization Unit	14
5 INSTALLING THE DETECTOR SYSTEM	16
5.1 Transport Considerations	16
5.2 Mounting	16
5.2.1 Mounting from Below	16
5.2.2 Mounting from Above	17
5.3 Grounding of the Detector	18
5.4 Connection to Dry Air or Nitrogen	19
5.5 Mounting the Detector Control Unit	19
6 TEMPERATURE AND HUMIDITY CONTROL	20
7 OPERATION PROCEDURE	21
7.1 Getting Started	21
7.2 Startup Procedure	21
7.3 Turning Off the Detector	21
7.4 Storing the Detector	22
7.5 Cleaning and Maintenance	22
8 TROUBLESHOOTING	23

DOCUMENT HISTORY

Current Document

Table 1: Current Version of this Document

Version	Date	Status	Prepared	Checked	Released
v1.6.1	2020-11-27	release	LW	SB, MM	LW

Changes

Table 2: Changes to this Document

Version	Date	Changes
v1.0.0	2017-04-09	First release.
v1.2.0	2017-09-04	EIGER2 R 500K integration.
v1.3.2	2017-09-04	PILATUS3 and EIGER2 R 500K API documentation integration.
v1.4.0	2018-06-19	EIGER2 Si detector series integration
v1.5.0	2020-03-02	New layout and rework of the technical specifications.
v1.5.1	2020-06-08	Fixed EIGER2 16M image of ground plate.
v1.6.0	2020-09-08	EIGER2 CdTe detector series integration.
v1.6.1	2020-11-27	EIGER2 S series integration.

1. GENERAL INFORMATION

1.1. Contact and Support





Address: DECTRIS Ltd.
Taefernweg 1
5405 Baden-Daettwil
Switzerland

Phone: +41 56 500 21 02
Fax: +41 56 500 21 01

Homepage: <http://www.dectris.com/>
Email: support@dectris.com

Should you have questions concerning the system or its use, please contact us via telephone, mail or fax.

1.2. Explanation of Symbols

Danger	#0
 Danger blocks are used to indicate immediate danger or risk to personnel or equipment.	
Warning	#0
 Warning blocks are used to indicate danger or risk to personnel or equipment.	
Caution	#0
 Caution blocks are used to indicate danger or risk to equipment.	
Information	#0
 Information blocks are used to highlight important information.	

1.3. Warranty Information

Caution

#1



Do not ship the system back before you receive the necessary transport and shipping information.

1.4. Disclaimer

DECTRIS has carefully compiled the contents of this manual according to the current state of knowledge. Damage and warranty claims arising from missing or incorrect data are excluded.

DECTRIS bears no responsibility or liability for damage of any kind, also for indirect or consequential damage resulting from the use of this system.

DECTRIS is the sole owner of all user rights related to the contents of the manual (in particular information, images or materials), unless otherwise indicated. Without the written permission of DECTRIS it is prohibited to integrate the protected contents in this publication into other programs or other websites or to use them by any other means.

DECTRIS reserves the right, at its own discretion and without liability or prior notice, to modify and/or discontinue this publication in whole or in part at any time, and is not obliged to update the contents of the manual.

2. USE OF THE EIGER2 S 9M

The EIGER2 S 9M detector system has been designed for the detection of X-rays produced by synchrotrons or laboratory sources. It is intended for indoor use only. For other applications, please contact DECTRIS technical support for additional information.

Caution

#2



Improper use of the DECTRIS detector system can compromise its safety and its functionality is no longer guaranteed.

Warning

#1



Do not use the detector in vacuum.

2.1. Product Return and Recycling

We recycle DECTRIS detector systems that are no longer suitable for use. If you are not using your DECTRIS detector system any more, send it back to us. We will make sure that your system is responsibly and safely recycled. This is free for customers who purchased a new DECTRIS detector system.

3. TECHNICAL SPECIFICATIONS

3.1. Specifications

3.1.1. Detector

Table 3.1: Technical Specifications

Number of modules (W x H)	3 x 6 = 18
Sensor	Reverse-biased silicon diode array
Sensor material	Silicon (Si)
Sensor thickness	450 μm
Pixel size (W x H)	75 μm x 75 μm = 5625 μm^2
Module size (W x H)	77.1 mm x 38.4 mm = 2961 mm ²
Pixel array format (W x H)	3108 pixel x 3262 pixel = 10 138 296 pixel
Active area (W x H)	244.65 mm x 233.1 mm = 57 027.915 mm ²
Intermodule gap	hor. 38 pixels, vert. 12 pixels
Image bit depth	32 bit
Readout bit depth	16 bit
Maximum count rate	1.7 $\times 10^9$ photons/s/mm ²
Adjustable threshold range	3.5 keV to 30 keV
Energy range	6 keV to 40 keV
Number of thresholds	two independent thresholds
Readout time	continuous readout, with zero dead time
Maximum frame rate ¹	40 Hz
ROI maximum frame rate	40 Hz
Point-spread function	1 pixel (FWHM)
Connection to control unit	4 x LC/UPC duplex connectors
Power supply	Internal power supply unit
Software interface	HTTP REST interface (via network connection)
Dimensions (W x H x D)	340 mm x 370 mm x 500 mm
Weight	42 kg
Overvoltage category	II
Pollution degree	II

¹ One frame per exposure, i.e. single-threshold or difference image

Table 3.1: Technical Specifications - continued

Maximum operating altitude	2000 m a.s.l.
----------------------------	---------------

3.2. Ratings

3.2.1. Detector

Table 3.2: Power Ratings

Detector power input	100 - 240 V AC, 50/60 Hz, 700 W C20 appliance inlet
<div>Warning #2</div> <div>⚠ Only use cables rated for 16 A.</div>	
Fuse	10 A slow-blow fuse. Type: Schurter 5x20 mm, 10 A, 250 V AC Part No. 0001.2514
<div>Warning #3</div> <div>⚠ Always replace fuses with the same type.</div>	
Detector external trigger input	High level: 2.1 – 5.0 V Low level: 0.0 – 0.8 V
<div>Caution #3</div> <div>⚠ Absolute maximum is 5 V. Applying a higher voltage will damage the detector.</div>	
External trigger input impedance	47 kΩ
Detector trigger output	High level: 2.3 V to 3.3 V Low level: 0.0 V to 0.6 V Max. current: 24 mA

3.2.2. Detector Control Unit

Information	#1
-------------	----



Please consult the user documentation of the DELL PowerEdge R940 for details.

Table 3.3: Detector Control Unit Ratings

Detector control unit power input	2 x 100V to 240V AC, 50/60 Hz, 5 A to 10 A, 750 W (Platinum) 1+1 redundant, hot swappable power supply unit
Dimensions (W x H x D)	482.4 mm x 130.3 mm x 776.46 mm
Weight	<50 kg
Chassis	3U

3.2.3. Thermal Stabilization Unit

Information	#2
-------------	----



Please consult the user documentation of the SMC HRS-012A thermal stabilization unit for details.

Caution	#4
---------	----



The maximum allowable coolant pressure is 3 bar.

Table 3.4: Thermal Stabilization Unit Ratings

	Value
Thermal stabilization unit power input	115 V version Single phase 100 VAC 50/60 Hz, 115 VAC 60 Hz, allowable voltage range $\pm 10\%$, 7.5 A (50 Hz) to 8.3 A (60 Hz) 230 V version Single phase 200 VAC to 230 VAC 50/60 Hz, allowable voltage range $\pm 10\%$, 4.6 A (50 Hz) to 5.1 A (60 Hz)
Dimensions (W x H x D)	377 mm x 615 mm x 500 mm
Weight	115 V version: 40.0 kg 230 V version: 43.0 kg
Typical flow	3 L min ⁻¹
Maximum operation pressure	3 bar

3.3. Ambient Conditions

Caution

#5



The EIGER2 S 9M detector is equipped with a temperature and a humidity sensor. When either sensor detects that the operating conditions are not met, the detector will shut off. However, as the sensors may not prevent damage, temperature and humidity should be monitored to avoid breaching the operation limits.

The EIGER2 S 9M detector is designed for indoor use only. The ambient conditions shown in table 3.5 must be satisfied. The stated values are for the ambient conditions.

Values inside the detector, in particular due to the dry-air or nitrogen supply, are different. These are described in section 5.4 and chapter 6.

Table 3.5: Detector Operating Ambient Conditions

Ambient Condition	Value
Operating temperature	+20 °C to +35 °C
Operating humidity	<80 % at 20 °C, non-condensing
Storage temperature	+15 °C to +40 °C
Storage humidity	<40 % at 20 °C, non-condensing

Caution

#6



Please consider the following points when storing the detector

- Make sure the temperature and the humidity inside the transport box does not exceed the specified range (use of a drying agent is required).
- Ensure that no condensation moisture develops if the detector is stored at low temperature.

4.1.2. Front Side of the Detector

Danger

#1



Danger of electric shock. Do not touch the Mylar® foil. The sensors behind the Mylar® foil are operated at high voltages. Touching the Mylar® foil can cause an electrical shock.

Warning

#4



Do not touch the Mylar® foil to avoid damage of the sensors.

Caution

#7



The cover may not protect the detector from a direct beam.

The detector comes with a protective cover (1.5 mm, Steel) for the front window, which should only be removed during operation. The sensors are behind a 12 µm thick Mylar® (PET) foil coated with aluminum to protect them from humidity, dust and from being touched.



Figure 4.2: The EIGER2 S 9M Detector with the Cover Removed (front view)

4.1.3. Back Side of the Detector

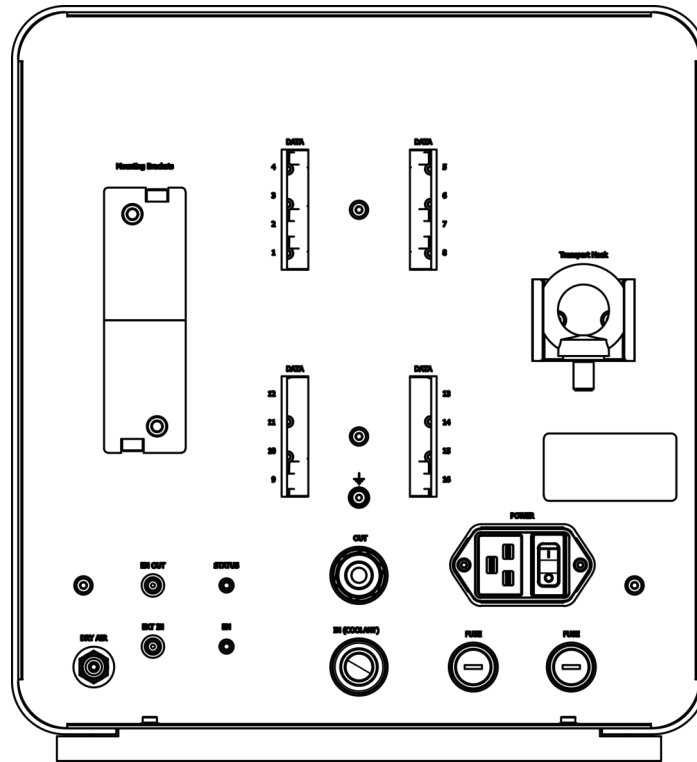


Figure 4.3: The EIGER2 S 9M Detector (back view)

4.1.4. Status LEDs

Table 4.1: The Meaning of the Status LEDs on the Detector Back Plane



LED	Behavior	Description
EN	Orange	Indicates the detector is in counting mode.
STATUS	Green steady	Detector running and hardware OK.
	Green blinking	Detector hardware OK, detector in Standby or Startup mode. ⇒ Press power button to turn on the detector.
	Red blinking	Detector may be overheating or humidity too high ⇒ Check cooling system and dry air supply
	Off	Detector has no power ⇒ Check the power adapter and mains connection.

4.1.5. Connectors and Connecting Cables/Pipes

Table 4.2: Electric Connectors and Connecting Cables

Connector	Description
DATA	<p>4 x LC/UPC duplex connectors DATA 1 -> detector control unit port 1 DATA 5 -> detector control unit port 5 DATA 9 -> detector control unit port 9 DATA 13 -> detector control unit port 13</p> <p>The unused data ports are closed with an EMI plug. Do not remove the plugs. Use Single Mode fiber optic patch cable with LC/UPC duplex connectors at both ends. We recommend to use the optic patch cables that are supplied with the detector system. Detector and detector control unit are equipped with 10GBASE-LR Single Mode SFP+ optical transceivers. Do not replace the optical receivers, as proper function of the detector system cannot be guaranteed otherwise. Please contact support@dectris.com if you need replacement transceivers.</p> <div style="background-color: yellow; padding: 5px; display: flex; justify-content: space-between;"> Caution #8 </div> <div style="display: flex; align-items: center; margin-top: 10px;"> <p>There must be a 4 x LC/UPC duplex connectors point-to-point connection between detector and detector control unit.</p> </div>
POWER	AC power connector (see table 3.2)
EXT IN	External trigger input (see table 3.2) Use a Lemo® Type 00 (NIM/CAMAC) cable.
EN OUT	Enable out, high when counting is enabled. Use a Lemo® Type 00 (NIM/CAMAC) cable.

Table 4.2: Electric Connectors and Connecting Cables - continued

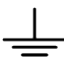Connector	Description
	Functional ground
<div>Information#4</div> <div> Although the detector might be already grounded via the mounting bolts, the detector should be grounded additionally via the functional ground connector at the back to establish a defined grounding.</div>	

4.2. Detector Control Unit

4.2.1. Configuration

Caution	#9
---------	----



Do not access or modify the operating system of the detector control unit.

The user interface of the detector control unit is accessible using a web browser. The detector control unit does not need any connections other than the power and Ethernet cables.

The detector control unit has to be connected point-to-point to the detector via 4 x LC/UPC duplex connectors. The detector control unit can be integrated into the site network infrastructure using one of the interfaces described in table 4.3. The detector control unit is optimized for performance and stability of operation. In order to achieve these goals we deliver the detector control unit with fixed firmware (BIOS etc.) and software (OS) version. The detector control unit must not be operated in an environment where unauthorized access is possible. The detector control unit does not provide authentication mechanisms and is not protected against malicious acts by unauthorized third parties.

Using the web front end, it is possible to restart the EIGER2 control service, trigger an update, and to shut down and to reboot the detector control unit. Any further control of the detector is carried out via the SIMPLON API (see separate documentation).



Figure 4.4: Front view of the detector control unit (symbolic picture).

Caution

#10



Pushing the power button on the front panel longer than 2 seconds will immediately halt the detector control unit. All image data on the detector control unit will be permanently lost.

Information

#5



Briefly pushing the power button on the front panel will shut down the detector control unit. May take up to 1 min.

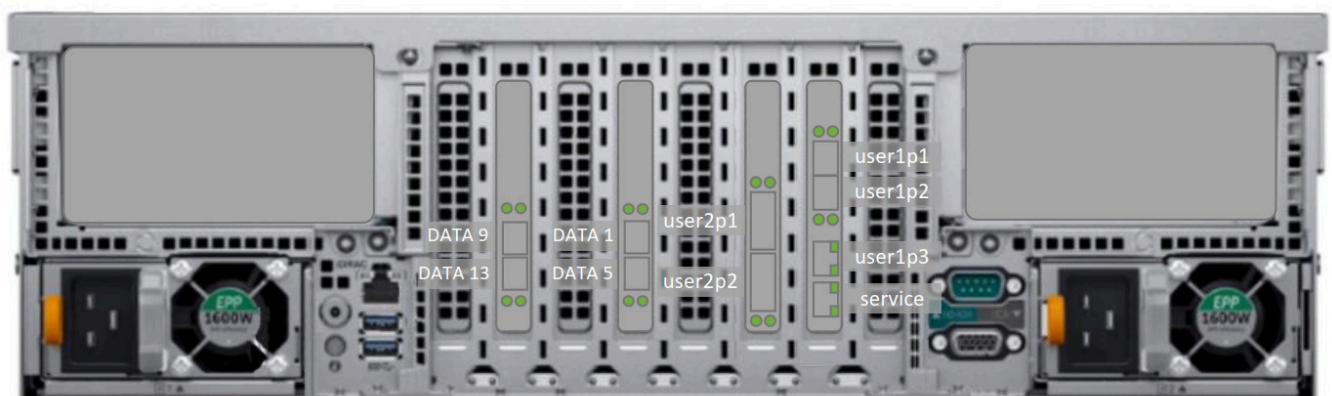


Figure 4.5: Back view of detector control unit with labeled network interfaces.

4.2.2. Connectors

Table 4.3: Detector Control Unit Connectors

Connector	Description
user1p1 (Integrated NIC)	Interface name: user1p1 (10 GbE SFP+) User configurable 10 GbE network interface Preconfiguration: DHCP
user1p2 (Integrated NIC)	Interface name: user1p2 (10 GbE SFP+) User configurable 10 GbE network interface Preconfiguration: Static 10.42.41.10 (Netmask 255.255.255.0)
user1p3 (Integrated NIC)	Interface name: user1p3 (1 GBase-T) User configurable GbE network interface Preconfiguration: DHCP
service (Integrated NIC)	Interface name: service (1 GBase-T) Fallback GbE network interface Preconfiguration: Static 169.254.254.1 (Netmask 255.255.255.0)
user2p1 (Slot 3)	Interface name: user2p1 User configurable 100 Gb QSFP28 network interface
user2p2 (Slot 3)	Interface name: user2p2 User configurable 100 Gb QSFP28 network interface
DATA	Detector interface ports DATA 1 -> detector control unit port 1 DATA 5 -> detector control unit port 5 DATA 9 -> detector control unit port 9 DATA 13 -> detector control unit port 13
2 x Power	AC Connector (redundant power supply)

See DELL owner's manual for further details.

4.3. Thermal Stabilization Unit

A thermal stabilization unit is required for the operation of the EIGER2 S 9M detector system.







The hoses and the detector are equipped with self-sealing valves to avoid dripping when connecting or disconnecting the tubes.

The tubing should be kept as short as possible to ensure the best flow.

Table 4.4: Operating Conditions

Condition	Definition
Operating temperature	The thermal stabilization unit has to be set to a temperature of 23 °C for normal operation.
Maximum operating pressure	3 bar

Table 4.4: Operating Conditions - continued

Condition	Definition
Coolant	Use 66 % distilled water and 34 % ethylene glycol.
	<div>Danger #2</div> <div>  Ethylene glycol can be seriously harmful to your health or fatal if handled incorrectly. Consider the packaging and safety instructions provided by your local supplier. </div>
<div>Information #6</div> <div>  Before operating the thermal stabilization unit, please read the User Manual of the thermal stabilization unit. </div>	
Caution	#11
	When connecting or disconnecting the cooling hoses, turn off the detector and the thermal stabilization unit.
Caution	#12
	When operating the detector, the thermal stabilization unit must always be turned on and the pump has to be activated (see user documentation of thermal stabilization unit).
Caution	#13
	Use opaque hoses to avoid the growth of algae.
Caution	#14
	Do not set the temperature of the thermal stabilization unit below the recommended operating temperature. Condensing moisture can develop and damage the detector.

5. INSTALLING THE DETECTOR SYSTEM

5.1. Transport Considerations

Warning

#5



Avoid vibration and shock when moving the detector.

Caution

#15



Use the included lifting eye for transporting the detector.

The detector has been delivered in a robust transport box. Please keep this transport box for transport or storage purpose.

5.2. Mounting

Caution

#16



Do not place the detector near heat sources or in a place subject to direct sunlight, excessive dust or mechanical shock.

Make sure that the detector has adequate ventilation.

- Do not cover any air intakes or outlets.
- Place the detector in a location with adequate air circulation.
- Make sure the detector has enough space for proper ventilation (minimum wall distance: 100 mm).
- Do not operate the detector in a closed environment.

The detector can be mounted in the ways which are described below.

5.2.1. Mounting from Below

Warning

#6



It is strictly forbidden to add any threads to the detector base plate or to the detector housing.

The detector should be mounted using the four internal M6x1 threads as shown in figure 5.1 (indicated with yellow circles).

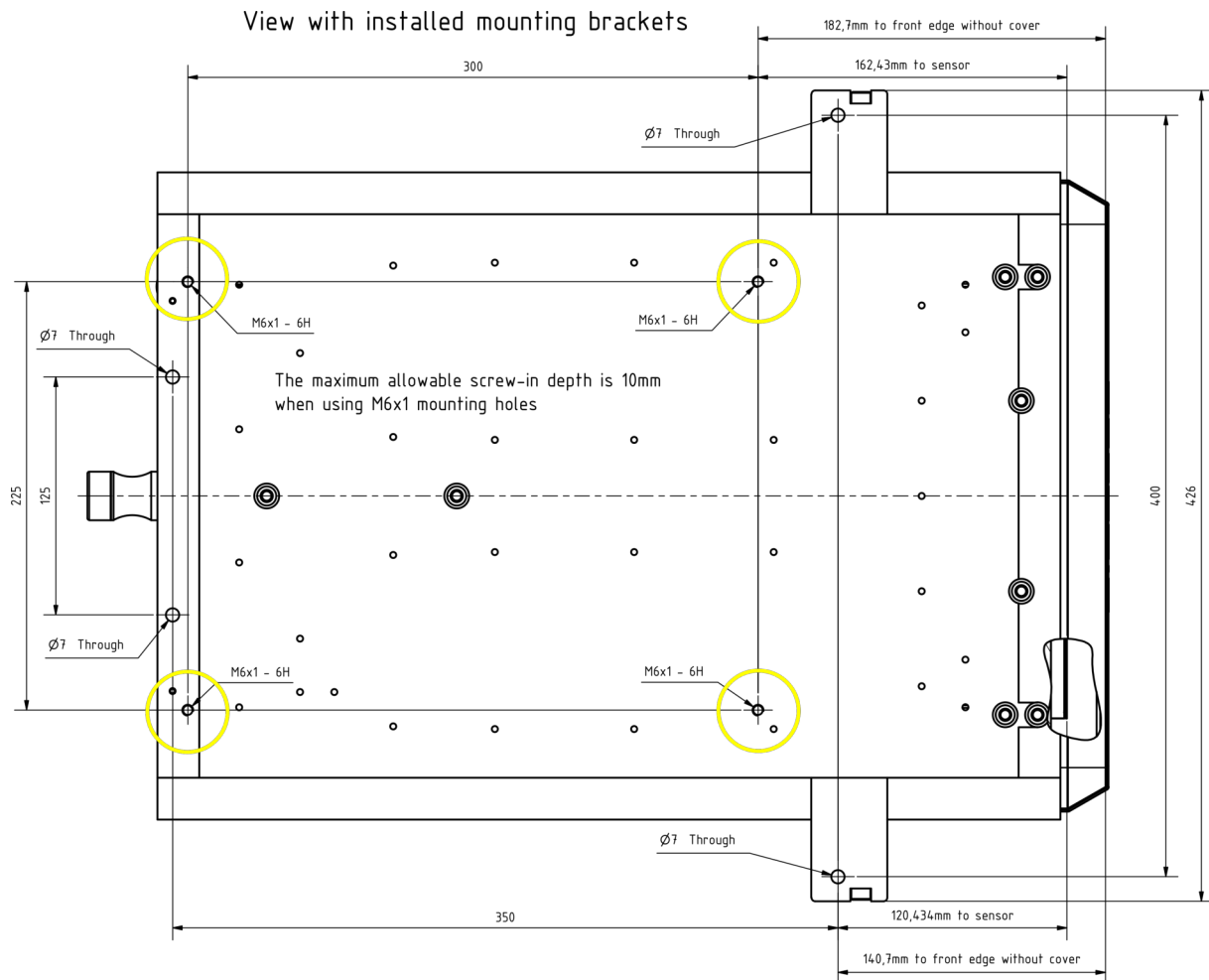


Figure 5.1: Drawing of the EIGER2 S 9M Detector Base Plate (also printed separately in the user documentation folder)

Caution

#17



The four M6 screws must not intrude into the detector more than 10 mm.

5.2.2. Mounting from Above

Use the mounting brackets as depicted in figure 5.2. These mounting brackets have to be mounted on the base plate of the detector. The detector should be mounted using all four outer 7 mm holes (indicated with yellow circles).

Caution

#18



Make sure the mounting brackets are mounted and properly tightened using the screws provided. The placement of the screws is indicated with diamonds in figure 5.2.

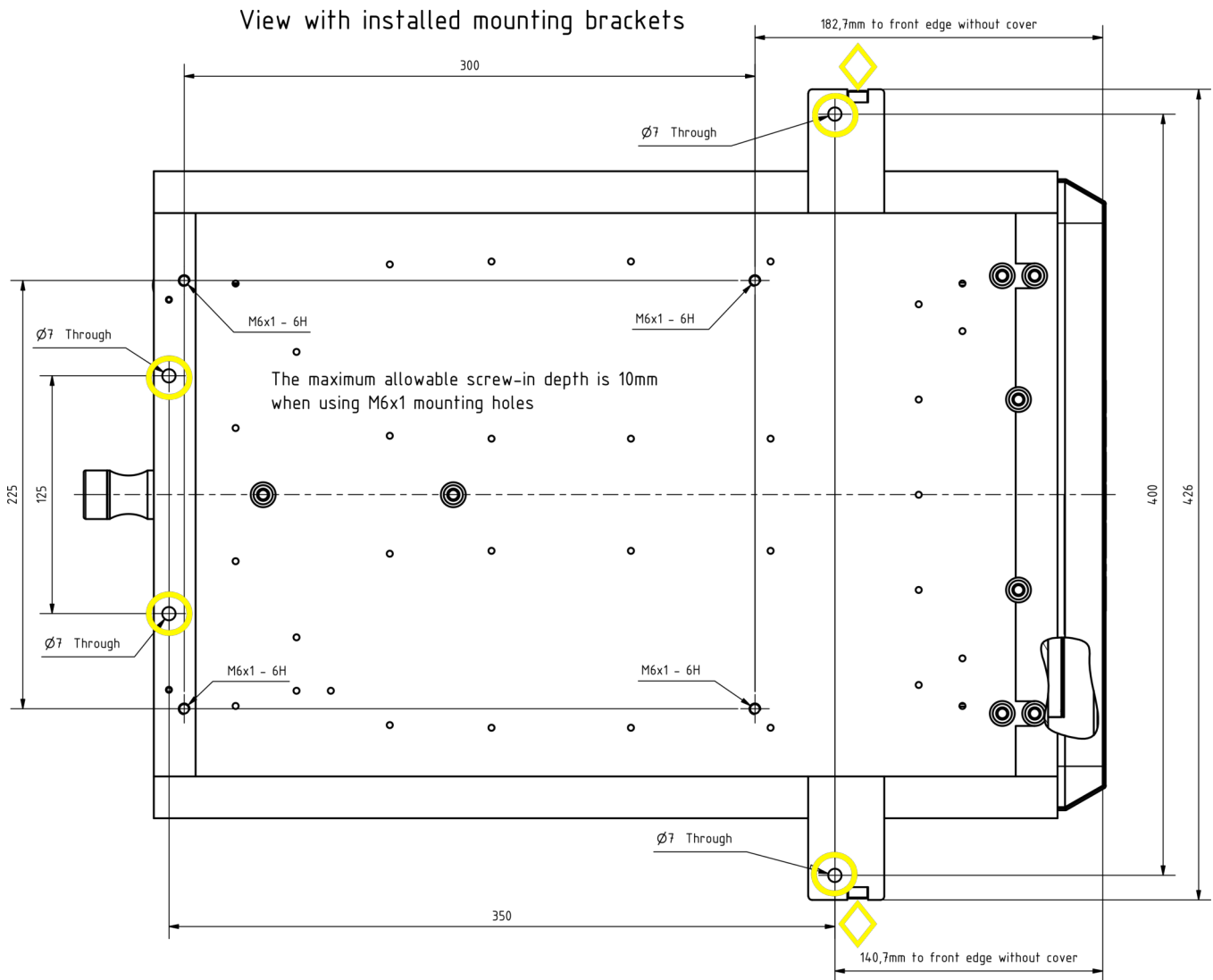


Figure 5.2: Drawing of the EIGER2 S 9M Detector Base Plate with Mounting Brackets (bottom view)

5.3. Grounding of the Detector

Caution

#19



The main plug of the detector control unit has to be connected to a grounded power outlet.

Although the detector might be already grounded via the mounting bolts, the detector should be grounded additionally via the functional ground connector at the top in order to establish a defined grounding. Although the detector might be already grounded via the mounting bolts, the detector should be grounded additionally via the functional ground connector at the back in order to establish a defined grounding.

5.4. Connection to Dry Air or Nitrogen

Warning

#7



When venting with nitrogen, take proper precaution against the risk of asphyxiation caused by oxygen displacement from nitrogen. Ensure sufficient ventilation and oxygen level monitoring. Use compressed air for venting large vessels or in confined spaces.

Caution

#20



Humidity might damage the detector. Make sure that the detector is operated within the allowed ambient conditions (see section 3.3).

The EIGER2 S 9M detector has to be connected to a dry air (or nitrogen) source to avoid humidity and condensation damage when it is outside of the storage box. For information on system connections, refer to the section 4.1.5 and for storage of the detector system refer to the section 7.4.

Information

#7



Oil free, dry air of <20 % relative humidity or nitrogen must be used.

- The recommended flow is 5 l h^{-1} to 10 l h^{-1} (at 2 bar).
- For reliable operation we recommend dry air of <5 % relative humidity.
- The gas pressure must not exceed 2 bar.
- The minimum gas pressure is 1 bar.
- The humidity control shuts down the power of the detector modules when the humidity is too high (see chapter 6).

5.5. Mounting the Detector Control Unit

Caution

#21



Make sure that the detector control unit has adequate ventilation.

The detector control unit can be mounted in a standard 19 inch rack, which has to be properly grounded.

6. TEMPERATURE AND HUMIDITY CONTROL

The EIGER2 S 9M detector has a combined temperature and humidity sensor.
The temperature and humidity control shuts down the detector when the humidity or the temperature of the sensor exceeds the following limits:

Table 6.1: Temperature and Humidity Limits

Shutdown Temperature	Shutdown Humidity
15 °C	35 °C
	40 % at operation / 40 % at start-up

The communication with the detector control unit will remain active after a temperature shut down (only power of the modules shuts down).

Warning

#8



The temperature and humidity control cannot prevent condensation issues and resulting damage to the sensor due to improper use. Always make sure that the detector is set to the correct temperature (thermal stabilization unit temperature set to 23 °C).

Information

#8



The detector has an internal thermal protection switch. If the thermal protection switch is activated, it turns the detector off to prevent damage. In this case all LEDs will be off and no communication is possible any more with the detector.

Information

#9



If the humidity is outside the specified range, the software will prevent operation.
The user can check the temperature and humidity via the API, as long as the temperature is not out of range. If the temperature breaches safe conditions the thermal protection switch will be triggered and the detector will completely switch off.

To start the detector correctly, please refer to section 7.1 and execute the correct startup procedure.

Make sure that the cooling unit is running at the recommended temperature (according to section 4.3) and that Nitrogen or dry air flow is turned on at the recommended flow rate, given in section 5.4. Then restart the software.

Information

#10



A free-flowing air stream is mandatory in order to properly cool the electronics inside the detector. Do not cover any ventilation holes.

7. OPERATION PROCEDURE

Before operating the detector, make sure you have read the Technical Specifications and the User Manual.

7.1. Getting Started

Before switching on:

- Mount the detector properly.
- Connect the detector to ground potential, using the functional ground connector.
- Connect the detector to power supply.
- Connect the detector to a nitrogen or dry air source, capable of supplying at least the minimum recommended flow rate.
- Connect the coolant hoses. Make sure they are properly mounted on both sides.
- Set the temperature to 23 °C on the thermal stabilization unit and turn it on. If the detector was not at room temperature, wait until the thermal stabilization unit has reached stable operation.
- Connect the power cable, the local network cable, and the detector data cable to the detector control unit. (If more than one data cable is required, please pay attention to the numbering of the cables as described in table 4.2)

7.2. Startup Procedure

Please use the following startup procedure:

- Turn on the dry air or nitrogen at least 30 min before turning on the detector.
- Turn ON the power switch at the back of the detector.
- Turn on the detector control unit. Wait at least 5 min before trying to connect.

The detector is now ready to use.

Information

#11



The software start-up procedure is described in detail in the User Manual.

7.3. Turning Off the Detector

To turn off the detector:

- Turn OFF the detector control unit.
- Do not remove the nitrogen/dry air connection. It is a requirement that it is left at the recommended flow rate according to section 5.4.

Warning

#9



As long as the detector power cable is connected, the detector has to be considered under power.

7.4. Storing the Detector

Information

#12



Even if the detector is not in operation, it is recommended that the dry air or nitrogen flow is maintained to reduce the risk of humidity damage to the detector.

Please follow these instructions:

- Put the detector in a plastic bag, add at least 200 g of drying agent (i.e. silica gel) into the bag and seal it air-tight.
- Check the humidity and change the drying agent frequently for compliance with the storage requirements in section 3.3.

7.5. Cleaning and Maintenance

Caution

#22



The Mylar® foil must not be touched or cleaned. If it is damaged, please contact DECTRIS technical support.

The housing can be cleaned with a soft tissue.

The EIGER2 S 9M detector does not require any maintenance.

Please refer to the user documentation of the thermal stabilization unit for detailed information about the maintenance of your thermal stabilization unit.

8. TROUBLESHOOTING

table 8.1 provides an overview of possible problems with the detector system and instructions in order to solve the problems. If the problem you are experiencing is not listed below or if the instructions do not help, please contact support@dectris.com.

The LEDs at the back of the detector can provide valuable information for troubleshooting. Check section 4.1.4 for further informations.

Table 8.1: Troubleshooting

Problem	Cause	Solution
Detector control unit does not start properly.	Detector control unit is not powered.	Check the User Documentation of the detector control unit (see section 3.2.2).
Communication error, the detector is not found at startup.	Data cable is not connected or defective.	<p>Check the connection between detector control unit and detector. Make sure that there is a direct, peer-to-peer connection between the detector control unit and the detector.</p> <p>Avoid tangling or strong bending of the Ethernet data cable.</p> <p>Check the status of the LINK LED. If the detector control unit and the detector are powered and correctly connected, the LINK LED should be green (Takes up to 30s after power up).</p>
Detector shuts down.	Temperature or humidity error.	<p>Check that the detector is properly supplied with coolant and check the temperature of the coolant at the front panel of the thermal stabilization unit.</p> <p>Check the flow of nitrogen or dry air.</p> <p>Check the temperature of the detector using the SIMPLON API and wait until the detector cools down.</p> <p>Restart the detector again.</p>
The detector fails to turn on.	The power cord is not connected or the plug is incompletely inserted.	Connect the power cord firmly. Check the LED on the external power supply.
	The fuse is blown.	Replace the fuse, see section 4.1.5.
	The temperature is over the critical limit. The thermal protection was triggered.	Check the thermal stabilization unit. The detector will power on again, as soon as the temperature is within the allowed operating conditions.
Image acquisition not possible.	Detector is not properly initialized.	Initialise the detector via the SIMPLON API. (See API Reference)

Table 8.1: Troubleshooting - continued

Problem	Cause	Solution
Detector housing is humid.	Ambient humidity around the detector exceeds the operating conditions.	Shut down the detector immediately and check the humidity. Power up the detector only when the ambient humidity has been reduced.
